

CETAC TECHNOLOGIES

SERVICE BULLETIN

Subject: Modification of ASX-500 home location flag.	Bulletin No:	SB0905
	Date Issued:	1/27/97
	Priority:	Low

Product: ASX-500

Serial Numbers: 019705ASX to current
Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)

Service Location: Customer Site CETAC Technologies

Responsibility: Customer Trained CETAC Representative

Replacement Parts: Customer CETAC Technologies

DESCRIPTION

Upon modification of the carriage assembly to improve the resistance to corrosive effects, the length of the home location flag has been increased by ¼ inch. This longer flag can not be used on autosamplers prior to serial numbers 019705ASX because the probe alignment will be incorrect. Autosamplers with serial numbers after 019705ASX can not use the shorter flag because the alignment will be incorrect.

IMPLEMENTATION

Each spare part kit that utilizes these flags will contain 3 of each length of the flags to insure that the correct flag is available. To determine the correct flag to use refer to the serial numbers as listed above. If you have any doubts concerning the correct length flag to use, contact your nearest CETAC service center.

PARTS

Part Number	Description
-------------	-------------

AVAILABLE DOCUMENTATION