

CETAC TECHNOLOGIES

SERVICE BULLETIN

| | | |
|--|---------------------|-----------|
| Subject: Autosampler X-Axis Sensor Board Replacement | Bulletin No: | SB1001 |
| | Date Issued: | 9/21/2004 |
| | Priority: | Medium |

Product: ASX-510HS, ASX-520

Serial Numbers: 470976 on X-axis board
Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)

Service Location: Customer Site CETAC Technologies

Responsibility: Customer Trained CETAC Representative

Replacement Parts: Customer CETAC Technologies

DESCRIPTION

Replacement of X-axis sensor board to upgrade 4-pin connector to 6-pin connector.

IMPLEMENTATION

Replace X-axis sensor by removing the two screws and putting the new sensor in its place.

The Y-axis sensor also needs to be replaced. See available documentation for instructions.

Connect the Y-axis sensor to the X-axis sensor.

PARTS

| Part Number | Description |
|-------------|---------------------|
| SP6107 | Y-axis sensor board |
| SP6106 | X-axis sensor board |

AVAILABLE DOCUMENTATION

SVC119.pdf Guide to Replacing the Y-Axis Sensor Board on the ASX-520 & ASX-520HS Autosamplers