

CETAC TECHNOLOGIES

SERVICE BULLETIN

Subject: Autosampler Main Board Replacement	Bulletin No:	SB1002
	Date Issued:	9/21/04
	Priority:	Medium

Product: ASX-510HS, ASX-520

Serial Numbers: Main board #BDC011R0
Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)

Service Location: Customer Site CETAC Technologies

Responsibility: Customer Trained CETAC Representative

Replacement Parts: Customer CETAC Technologies

DESCRIPTION

Replacement of the main board on the autosampler may cause the main board to malfunction if the X-axis sensor board is not replaced.

IMPLEMENTATION

To replace the main board, the X-axis sensor board needs to be replaced at the same time. See available documentation for instructions.

PARTS

Part Number	Description
SP6111	Main board

AVAILABLE DOCUMENTATION

SVC118.pdf Guide to Replacing the Main Board on the ASX-520 & ASX-520HS Autosamplers

SB1001.pdf Autosampler X-Axis Sensor Board Replacement