

# CETAC TECHNOLOGIES

## SERVICE BULLETIN

<b>Subject:</b> Autosampler Power Connector Defective	<b>Bulletin No:</b>	SB1003
	<b>Date Issued:</b>	9/21/04
	<b>Priority:</b>	Low

**Product:** ASX-260, ASX-410, ASX-510HS, ASX-520  
**Serial Numbers:** Prior to 100401A260, 100401A410, 100401A520  
*Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)*

**Service Location:**  Customer Site  CETAC Technologies

**Responsibility:**  Customer  Trained CETAC Representative

**Replacement Parts:**  Customer  CETAC Technologies

### DESCRIPTION

Autosampler power connector has an inner connector that is weak that can break in shipping or when pressure is applied to it.

### IMPLEMENTATION

Replace connector by removing screws and cutting the wires. Then solder wires to the new connector and putting the screws back on.

### PARTS

Part Number	Description
SP6104	Switch/power input kit

### AVAILABLE DOCUMENTATION