

CETAC TECHNOLOGIES

SERVICE BULLETIN

Subject: Guide shaft corrosion	Bulletin No:	SB1007
	Date Issued:	1/25/05
	Priority:	Medium

Product: ASX-510, ASX-520, ASX-520HS, ASX-260, ASX-130

Serial Numbers: ALL
Serial Numbers start by month (MM), year (YY) and units built that month (i.e., MMYyxx)

Service Location: Customer Site CETAC Technologies

Responsibility: Customer Trained CETAC Representative

Replacement Parts: Customer CETAC Technologies

DESCRIPTION

Corrosion of the guide shaft has been observed in acidic environments.

IMPLEMENTATION

Loosen set screws in shaft and rotate 5-10 degrees to give a new bearing surface. Tighten set screws.

Clean guide shaft with acetone. Apply thin layer of PTFE lubricant to guide shaft.

PARTS

Part Number	Description
SP7033	Super Lube, PTFE Based Lubricant

AVAILABLE DOCUMENTATION